



Software Version:
1.7.1.3055

Release Date:
6.23.2021

Contents

- Changes Since 1.7.0.2459 Release
 - Features
 - Updates
 - Known Issues
 - Installation instructions

1 Changes Since Last Release

1.1 Features

- Ability to hide ROI label in Array View legends
- Read-only rights for .sprdata file opening to allow in process runs to be viewed

1.2 Updates

- Peptide mapping analyte dendrogram incorrectly included ligand data
- Normalization bar in Epitope Mapping would display off-scale
- Peptide Mapping Heat Map sensorgrams data not updating to included/excluded options
- Epitope Binning bins being assigned incorrectly after sample name change.
- UI issue: no limits on the input fields (Binning, Analysis tabs)
- Incorrect icon of full-screen mode on the serial and Overlay views (Active+Ref, Ref Subtracted, Ref Only)
- Incorrect checkbox behavior in the Ligands table on the Data tab so selected/deselected data
- Data are updating incorrectly on the overlay view.
- Peptide Mapping and Mutant Mapping modes incorrect checkbox behavior on Data tab
- Incorrect behavior of the saving and reopening options
- Control antigen disappears while sorting by name (Inj ID) in the Analytes table for Mutant Mapping
- Error with Batch analysis in th Premix Binning
- Unstable state error appears after clicking paste data to main Analyte table in Peptide Mapping





- Heat map in Peptide and Mutant Mapping modules not included in export
- Error in sorting in the Analytes table Peptide and Mutant Mapping
- Heat map not exporting in analyzed Epitope Binning file
- When the lock normalization bars option is active, zoom on sensorgrams and curve selection is disabled

1.3 Known Issues

- None

2 Installation/Upgrade Instructions

- i. Run the Epitope installer executable file
- ii. Accept User Account Control permissions
- iii. Complete standard installation steps as instructed
- iv. Open application from Windows menu or desktop shortcut

