



Software Version:  
1.8.0.3603

Release Date:  
06.17.2022

## Contents

- Changes Since 1.7.1.3055 Release
  - New Features
  - Updates
  - Known Issues
  - Installation instructions

### 1 Changes Since Last Release

#### 1.1 New Features

- Improved performance for file opening, saving, and exporting
- Improved performance on Data tab
  - Link Group reassignment
  - Column sorting
  - General navigation and actions
- Improved error reporting interface

#### 1.2 Updates

- Array View tab empty if one ligand or analyte group is selected
- Assigning to Complex and assigning to Excluded options aren't working in Steady State > array view page
- Tables are empty in Kinetic Snapshot and Quant tabs in the exported file
- Data isn't dynamically updated on the Kinetics > Array View page
- Link Group reassignment issue: Serial View/Overlay View charts not fully displayed after reassignment
- Incorrect copying of data from the Ligands table in the LSA files
- Chart legend isn't displayed in Report Point > Quant
- The chart is shifted along the X-axis if the linear fitting algorithm is applied several times on Quant tab
- The application crashes after checking the "Display Units" checkbox on Quant
- IsoAffinity Values in Export file all have default values
- The Kinetic Analysis table is exported incorrectly when the "Show error ()" checkbox is unchecked (kd ka km mode)





- The exported file has extra data on the "Kinetic Snapshot" tab
- Export error appears after applied kd ka km fitting in Kinetics > Kinetic Analysis
- Leading blank option not functioning correctly
- Incorrect behavior of excluding ligands on the Analysis tab
- Report Point not working for Surface Prep Lawn
- Default groups don't appear in the list on the Analysis tab > Stats subtab
- After removing ligand sets and opening a new file, an error appears
- Errors when exporting to Excel
- Chart from Steady State > Array View does not display in exported file
- Array View tab empty if one ligand or analyte group is selected
- Tables from Quant export do not match those in the application
- Standards outside of Min/Max allowable range are excluded from fitting and plotting
- Report Point -> Quant R-squared Value uses the standard values not experimental values
- An unstable error appears if manually updating standard concentration after saving and reopening the file
- Error after manually updating standard concentration values used in Fitting on Quant Page
- .kitx file opening with no concentration values for Unknowns in Quant
- Array View of exported steady state data is missing sensorgrams
- An unstable error appears after using "Copy Data To Clipboard" option (Analyte data)

### 1.3 Known Issues

- Flagged Data isn't included in Analysis with the "Use Nickname" option
- Chart Text Size doesn't change correctly on Steady State > Array View
- Kinetics tab: Copy to Clipboard Function is Missing X-Axis Labels and Title

## 2 Installation/Upgrade Instructions

- i. Run the Kinetics installer executable file
- ii. Accept User Account Control permissions
- iii. Complete standard installation steps as instructed
- iv. Open application from Windows menu or desktop shortcut

