



# Increasing DNA Encoded Library (DEL) Screening Resolution Using HT-SPR

## Key Takeaways

- Detailed kinetics for >5,000 compounds per week
- Compatible with DEL formatted fragments, small molecules, TPDs, macrocycles, etc.
- Increase confidence in hits prior to off-DNA scale up

## Introduction

DNA encoded library (DEL) technology has permitted substantial leaps in compound screening by enabling a more facile way to interrogate libraries and develop lead compounds<sup>1</sup>. While a huge benefit to DEL technology is its reduction in the sheer number and tracking of discrete compounds through the screening process, it does come with a limitation in binding properties that can be gleaned for potential hits. In practice, from an initial screen numbering millions, the resulting thousands of hits are then reduced to around 50 compounds carried forward into downstream assays with limited characterization of their binding properties<sup>2,3</sup>. These fundamental properties include binding kinetics as well as target epitope. To address this resolution gap in DEL screening, demonstrated here is a technique to further characterize hits with improved resolution using high-throughput surface plasmon resonance (HT-SPR). HT-SPR characterizes the full kinetic profile for both weak and strong binders. Assay resolution can be further

increased by inclusion of high-homology family members, truncated or mutated forms, binding partner disruption, etc., yielding a richer profile of lead candidates. With a screening capacity of thousands per week, HT-SPR affords a high-resolution technique that matches

the throughput needs in this phase of discovery between the full library and a handful of leads. Additionally, this approach can work with any moiety having DNA attachment, including screening for targeted protein degraders (TPDs) and macrocycles.



**Figure 1: The Carterra LSA**

## Kinetic Screening Thousands of DEL Compounds

The Carterra LSA (**Figure 1**) is a unique technology for HT-SPR screening of binding interactions in

real time. Novel microfluidics allow up to 384 samples to be arrayed on the biosensor surface, followed by screening in parallel against hundreds of injected samples. Up to 150,000 interactions can be assessed in a single run depending on the assay design. Data analysis is highly flexible using either dedicated software or else by interfacing with third-party data analysis packages.

**Figure 2** highlights the basic steps involved in creating an array of up to 384 DEL compounds. Key to enabling this approach is the use of a DNA tag that encompasses a region capable of hybridizing with the corresponding locked nucleic acid (LNA) surface. The same DNA tag could also include a coding region. The surface can easily be stripped of the DELs using a simple mixture of salt and base, and a new array then captured. The LSA instrument has plate capacity for up to 1,152 DEL compounds in a single run and allows for comprehensive kinetic screening of >5,000 compounds per week.



**Figure 2:** HT-SPR workflow for high resolution measurement of DEL kinetic affinities

### Kinetic Screening Results

Representative data for a detailed kinetics screen of 96 DEL compounds captured at four densities binding a 25 kDa protein target are shown in **Figure 3**. Colored highlighting of select interaction tiles is done by automatic data flagging features in the Kinetics analysis software which indicate low binding responses or deviations from a 1:1 Langmuir model. This approach not only stratifies binders versus non-binders, but also gives highly reliable measures of association and dissociation kinetics for thousands of compounds. **Figure 4** highlights a particular compound captured at 400, 200, 100, and 50nM onto the LNA chip surface. By having the capacity to capture at multiple densities in a single run, the effect of density on observe binding kinetics can be understood.



**Figure 3:** Tile view of sensorgram binding profiles for individual DELs. Red lines are 1:1 Langmuir model. Poor or complex binders are highlighted in gray or purple, respectively.

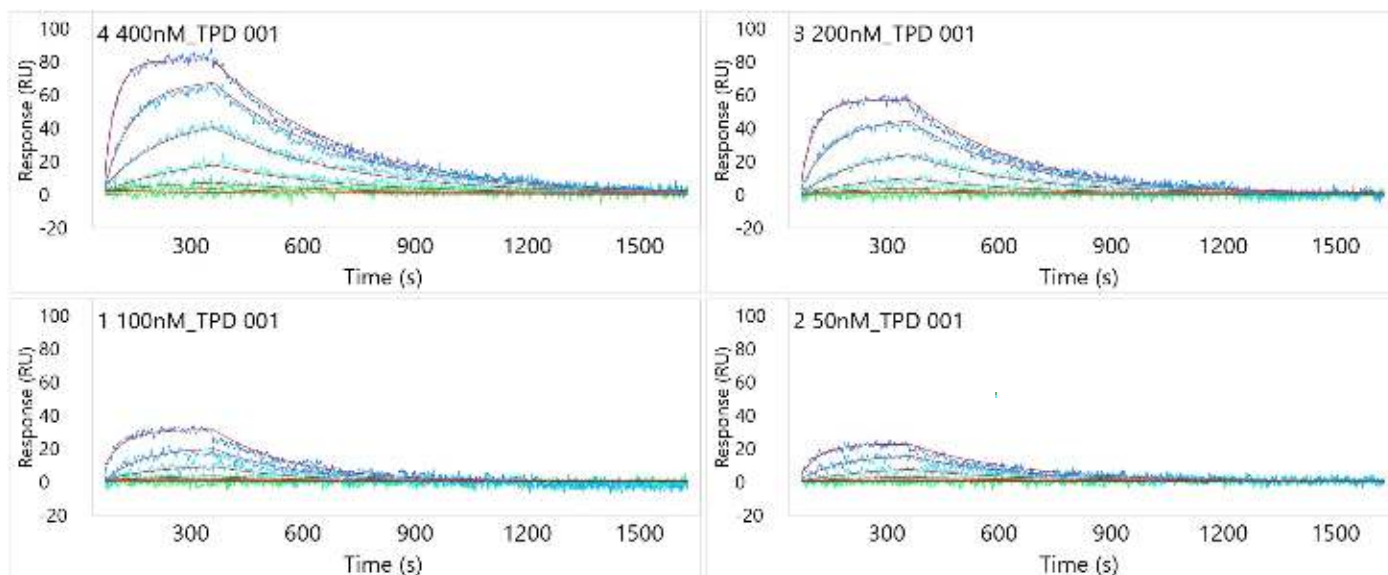


Figure 4: Kinetic profiles for a single DEL compound captured at four densities against a 25 kDa protein target.

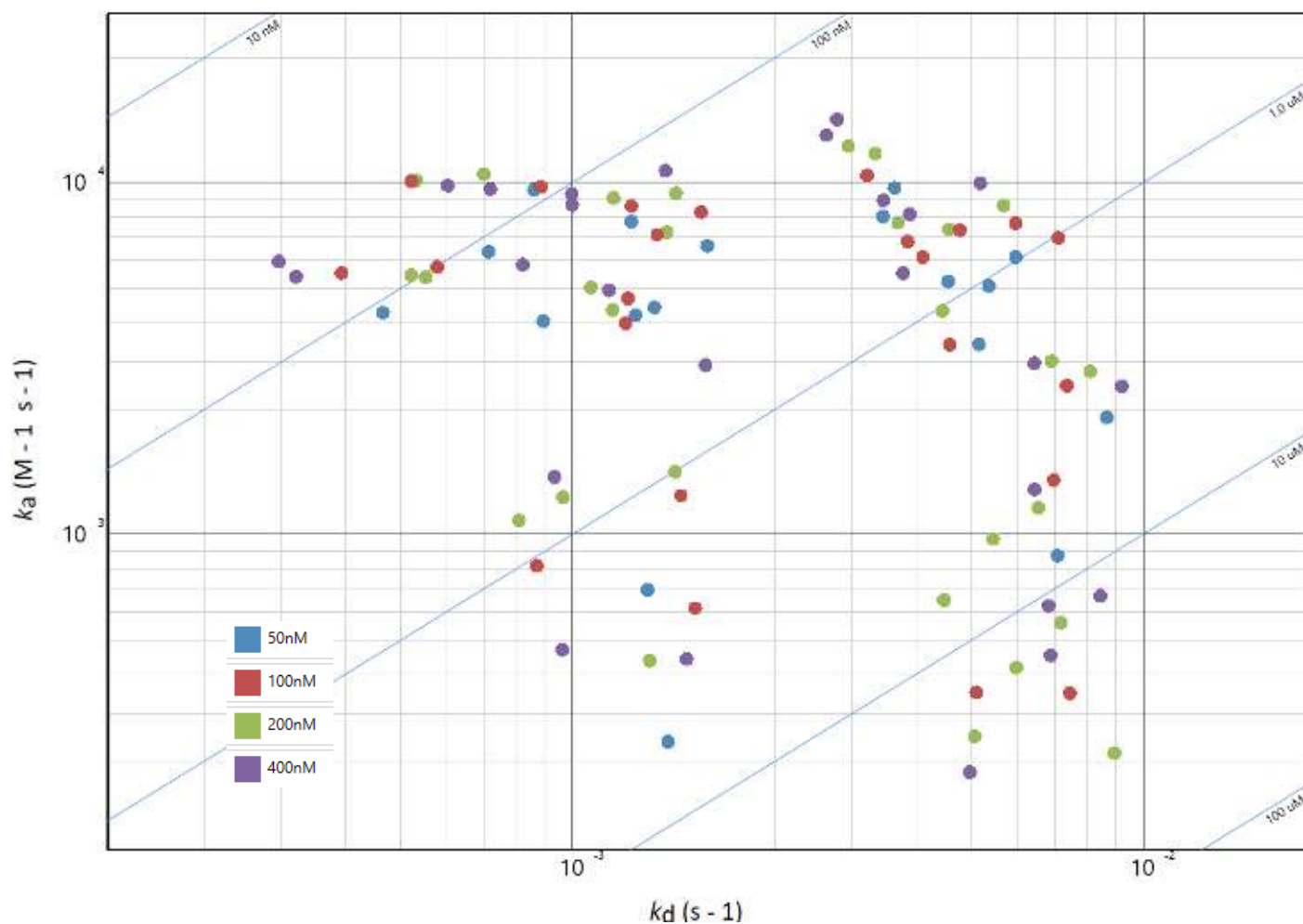


Figure 5: Isoaffinity plot of DEL compounds showing distribution of kinetics and colored by capture density.

## Highlights

Demonstrated here is an approach to DEL compound screening that seeks to bridge the gap between initial hit screening and off-DNA lead compound scale up. By increasing the data resolution prior to off-DNA screening, this approach affords a better chance of finding drug candidates with optimal properties from thousands of potential hits. While fragments and small molecules are typically the formats screened using DEL technology, this same assay can be applied to any molecule attached to DNA, including TPDs and macrocycles. Additionally, the assay has versatility beyond kinetic analysis where, for example, target isoforms and off-target reagents can be readily screened. Competitive assay formats are also adaptable to this approach and allow additional mechanistic questions to be addressed.

## References

1. Shi B, Zhou Y, Li X. Recent advances in DNA-encoded dynamic libraries. *RSC Chem Biol*. 2022 Feb 17;3(4):407-419.
2. Plais L, Scheuermann J. Macrocyclic DNA-encoded chemical libraries: a historical perspective. *RSC Chem Biol*. 2021 Oct 29;3(1):7-17.
3. Gironda-Martínez A, Donckele EJ, Samain F, Neri D. DNA-Encoded Chemical Libraries: A Comprehensive Review with Successful Stories and Future Challenges. *ACS Pharmacol Transl Sci*. 2021 Jun 14;4(4):1265-1279.

Carterra technology is protected by the following patents and other patents pending:  
8,210,119, 8,211,382, 8,383,059, 8,999,726, 9,682,372, 9,682,396, 10,825,548

© 2022 Carterra Inc. Carterra, the Carterra Logo and all other trademarks are property of Carterra unless otherwise specified.

AN114-REV006



info@carterra-bio.com

Headquarters:  
825 N 300 W Suite C309  
Salt Lake City, UT 84103

[www.carterra-bio.com](http://www.carterra-bio.com)